



CHECKLIST FOR WHEN YOU ARE DONE SCANNING

<p style="text-align: center;">Have you logged into the blue book in the control room?</p> <ul style="list-style-type: none"> Record date& time of scanning, PI's name, project title, user name
<p style="text-align: center;">Have you turned off the anesthesia vaporizer?</p> <ul style="list-style-type: none"> Applies to the MR and animal preparation room!
<p style="text-align: center;">Have you disconnect all anesthesia gases?</p> <ul style="list-style-type: none"> Disconnect the white O₂ and blue N₂O lines from the wall. If left unplugged, the user will be charged for the leaked gas! Applies to the MR and animal preparation room!
<p style="text-align: center;">Have you turned off the hot water circulation system?</p> <ul style="list-style-type: none"> If removing the bed base, disconnect the water lines from the base
<p style="text-align: center;">Have you disconnected the battery from the temperature module and connected it to the charging port on the respiration module?</p> <ul style="list-style-type: none"> The batteries deteriorate if left half-charged
<p style="text-align: center;">Have you disconnected the RF coil?</p> <ul style="list-style-type: none"> Disconnect the RF coil from back of the magnet and place it on the shelf Handle all equipment gently and with care!
<p style="text-align: center;">Have you turned off the power in the cabinet with high power electronics?</p> <ul style="list-style-type: none"> FIRST, press the top (on/off) button, under <i>local</i> and above <i>h.v.</i> ONCE ALL THE LIGHT ON THE CABINET BELOW TURN OFF, switch off the system by pressing the green switch labeled <i>power</i>
<p style="text-align: center;">Have you cleaned the animal bed and the table?</p> <ul style="list-style-type: none"> Spray with liquid in the red bottle, leave acting for 10 min, wipe off Spray with water in the blue bottle and wipe off Think of the user scanning after you and his/her animals!
<p style="text-align: center;">Have you removed and stored all the chemicals (contrast agent, drugs)?</p> <ul style="list-style-type: none"> Store chemicals (except isoflurane!) and biological samples in the fridge (the lock combination is 451)
<p style="text-align: center;">Have you tidied up behind you?</p> <ul style="list-style-type: none"> Remove all syringes, catheters, contrast agents, etc... Use the group cabinet(s) in the storage room next to the animal preparation room
<p style="text-align: center;">Have you switched off the light in the MR room?</p> <ul style="list-style-type: none"> Light source in the MR room heats up the room
<p style="text-align: center;">Have you locked the doors behind you?</p> <ul style="list-style-type: none"> Lock the main door the MR facility